

Sartobran P 0.45 µm

Bioburden and Particle Reductive Filter Cartridges



Description

Sartobran P 0.45 µm rated filter cartridges are ideally suited for bioburden and particle removal from bio-pharmaceutical solutions for protection of subsequent downstream processing equipment or sterilizing grade filters. The unique low unspecific binding capacity of the cellulose acetate membranes assures highest protein yield and rapid preservative recovery.

Applications

Sartobran P filters are ideally suited for prefiltration of high value biological solutions and pharmaceuticals sensitive to adsorption as well as for final filtration of LVP's and Buffers. Typical applications are filtration of:

- Coagulation Factors, Albumine, IgG
- Bacterial and Viral Vaccines
- MAB's
- Bio-processed Pharmaceuticals
- Diagnostics
- Purified Protein Solutions
- LV P
- Buffers

Highest Product Yield

The cellulose acetate membrane of the Sartobran P filters provides the lowest unspecific adsorption of all membrane materials available for highest product recovery rates.

Performance

Due to the "built-in prefiltration" by a 0.65 µm membrane, Sartobran P 0.45 µm filters provide excellent total throughputs and higher flow rates at low differential pressure for gentle product treatment.

Flexibility

Sartobran P 0.45 µm filters are available in traditional cartridge formats and disposable capsules from 150 cm² to 1.8 m² for simple linear scale up and process flexibility.

Quality Control

Each individual element is tested for integrity by diffusion and bubble point test prior to be released assuring absolute reliability.

Documentation

Sartobran P cartridges are designed, developed and manufactured in accordance with an ISO 9001 certified Quality Management System. A Validation Guide and Extractables Guide are available for compliance with regulatory requirements.

Specifications

Materials

Prefilter Membrane:	Cellulose Acetate
Endfilter Membrane:	Cellulose Acetate
Support Fleece:	Polypropylene
Core:	Polypropylene
End Caps:	Polypropylene
O-Rings:	Silicone (optional EPDM or Viton)

Pore Size

0.65 µm + 0.45 µm

Available Sizes | Filtration Area

Cartridges | MaxiCaps

Size 1	10"	0.6 m ²	6.5 ft ²
Size 2	20"	1.2 m ²	12.9 ft ²
Size 3	30"	1.8 m ²	19.4 ft ²

Capsules | Mini Cartridges

Size 5	0.03 m ²	0.3 ft ²
Size 7	0.05 m ²	0.5 ft ²
Size 8	0.1 m ²	1.1 ft ²
Size 9	0.2 m ²	2.2 ft ²
Size 0	0.45 m ²	4.8 ft ²

Available Adapters Cartridges

21, 25, 27, 28

Available Adapter Mini Cartridges

15

Available Connectors Capsules | MaxiCaps

SS, SO, OO

Operating Parameters

Max. Allowable Differential Pressure:	5 bar 75 psi at 20 °C (Cartridges)
	4 bar 58 psi at 20 °C (Capsules)
	3 bar 43.5 psi at 20 °C (MaxiCaps)
	2 bar 29 psi at 80 °C (Cartridges and Capsules)
Max. Allowable Back Pressure:	2 bar 29 psi at 20 °C

Membrane Filter Cartridges

Specifications

Extractables

Sartobran P 0.45 µm rated filter cartridges meet, or exceed the requirements for WFI quality standards set by the current USP.

Regulatory Compliance

100% Individually integrity tested

Integrity test correlated to HIMA/ASTM F 838-83 Bacteria Challenge Test

Non-pyrogenic according to USP Bacterial Endotoxins

Passes USP Plastics Class VI Test

Non-fiber releasing according to 21 CFR

Sterilization

In-Line Steam Sterilization

134 °C, 20 min. at max differential pressure of 0.5 bar | 7 psi

Note:

Capsules and MaxiCaps cannot be in-line steam sterilized.

Autoclaving

134 °C, 2 bar | 29 psi, 30 min

Sterilization Cycles

In-Line Sterilization: Min. 25

Autoclaving: Min. 25

Technical References

Validation Guide:

SPK 5726-e


Extractables Guide:

SPK5720-e

Ordering Information


Order Code	Pore size [µm]
Standard Cartridges	
523**06D1----P	0.45
523**06D2----P	0.45
523**06D3----P	0.45
Capsules	
5231306D5--00--B	0.45
5231306D7--**--B	0.45
5231306D8--**--B	0.45
5231306D9--**--A	0.45
5231306D0--**--V	0.45

Water Flow Rates for Standard Cartridges and MaxiCaps

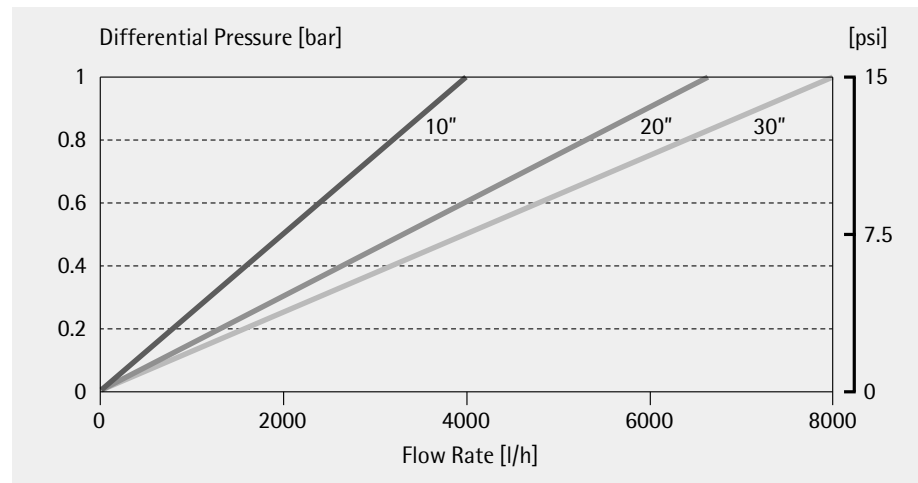

Standardized at 20 °C

Total Throughput Comparison

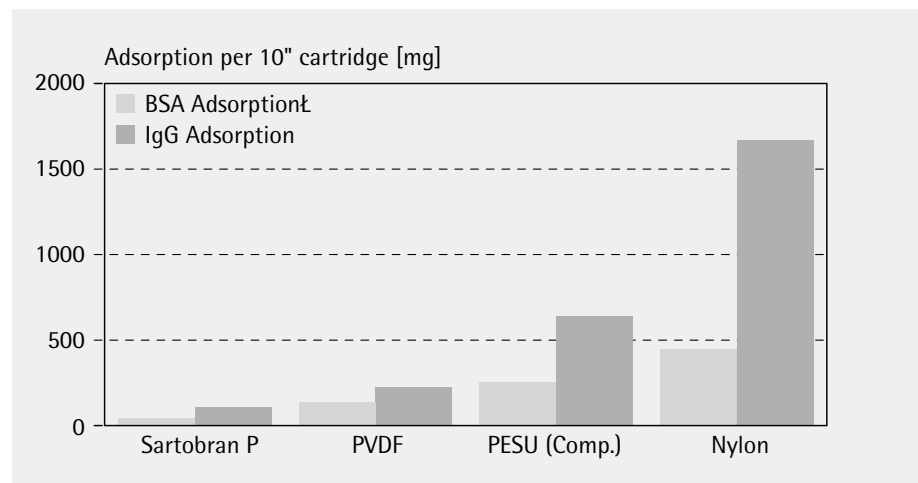

10" Cartridge format

Order Code	Pore size [µm]
MaxiCaps	
5231306D1--**	0.45
5231306D2--**	0.45
5231306D3--**	0.45
Mini-Cartridges	
5231506D7-----B	0.45
5231506D8-----B	0.45
5231506D9-----B	0.45

Sartorius AG
Weender Landstrasse 94-108
37075 Goettingen, Germany

Phone +49.551.308.0
Fax +49.551.308.3289
www.sartorius.com

USA +1.631.2544249
UK +44.1372.737100
France +33.1.69192100
Italy +39.055.634041
Spain +34.91.3586100
Japan +81.3.33293366

Specifications subject to change without notice.
Printed and copyrighted by Sartorius · W/sart-000 · G
Publication No.: SPK2050-e02121
Order No.: 85030-516-92